


PathSuite™ Professional Quick Start Guide

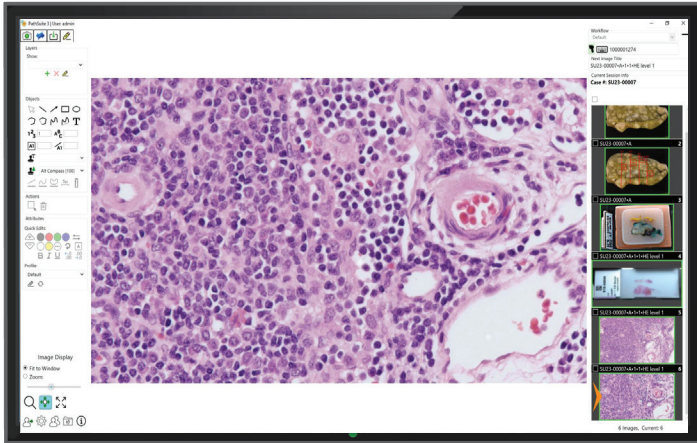


866-604-SPOT • support@spotimaging.com
spotimaging.com


PathSuite™ Professional Quick Start Guide



NOTE: Before beginning the installation process ensure that Windows is not anticipating a restart (*such as a Windows Updates*). The computer must be connected to the internet. If these requirements are met, double click the **PathSuite Professional Installer Icon**.

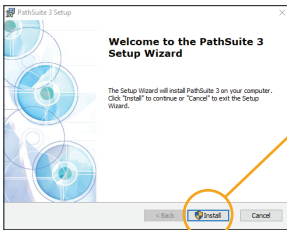


① Click the **“Install”** button and agree to any prompts

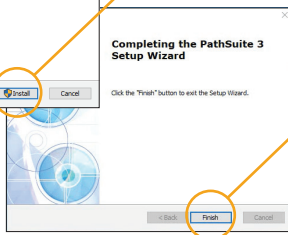


② Click the **“Finish”** button to exit the Installer.
NOTE: a computer restart is required!



③ Launch the **“License Manager”** via the desktop shortcut

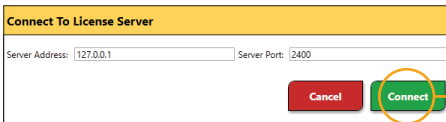


④ Confirm the Server Address is set to **“127.0.0.1”** and the Server Port is set to **“2400”**
Click the **“Connect”** button

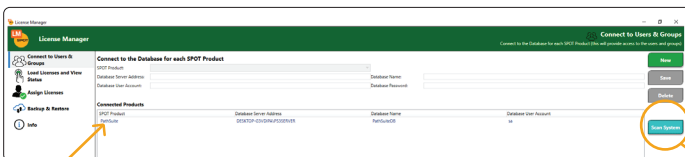


⑤ In the **“Connect to Users & Groups”** panel, press the **“Scan System”** button and confirm **“PathSuite”** is listed under **“Connected Products”**



- 6 In the **"Load Licenses and View Status"** panel, click the **"Load"** button and select your .lic file

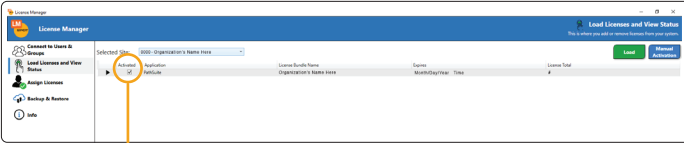

- 7 Wait for the license file to load and to automatically be activated (noted by a checked box in the **"Activated"** column)

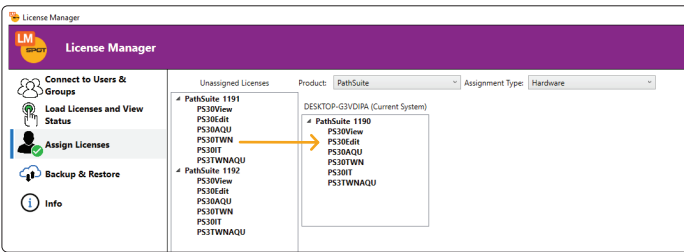

- 8 In the **"Assign Licenses"** panel, drag and drop a License from the **"Unassigned Licenses"** list to the **"Current System"**

- 9 Close License Manager



- 10 Launch **PathSuite** using the desktop shortcut

- 11 Log in for the first time with the following credentials,
Username: **admin** • Password: **admin**

- 12 During the first login the user will be prompted for the License Server Information. Confirm the Server Address is set to **"127.0.0.1"** and the Server Port is set to **"2400"**. Click the **"Save"** button

- 13 It is recommended to change the admin's account password using the **"Users"** button in the bottom left of the **PathSuite** application.