




10 Launch **PathSuite** using the desktop shortcut

11 Log in for the first time with the following credentials,  
Username: **admin** • Password: **admin**

12 During the first login the user will be prompted for the License Server Information. Confirm the Server Address is set to **"127.0.0.1"** and the Server Port is set to **"2400"**. Click the **"Save"** button

13 It is recommended to change the admin's account password using the **"Users"** button in the bottom left of the **PathSuite** application.


## PathSuite™ Professional Quick Start Guide

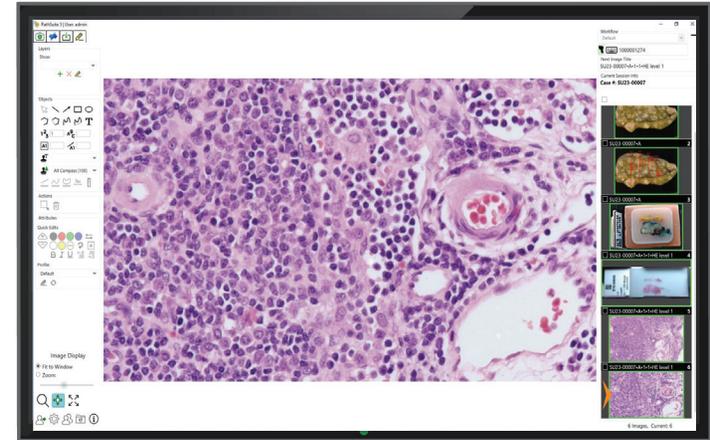


# PathSuite™ Professional Quick Start Guide



**NOTE:** Before beginning the installation process ensure that Windows is not anticipating a restart (such as a Windows Updates). The computer must be connected to the internet. If these requirements are met, double click the **PathSuite** Professional Installer Icon.

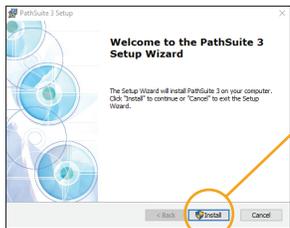


1 Click the **"Install"** button and agree to any prompts

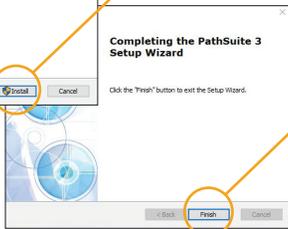


2 Click the **"Finish"** button to exit the Installer.  
**NOTE:** a computer restart is required!




3 Launch the **"License Manager"** via the desktop shortcut

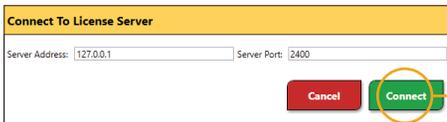

4 Confirm the Server Address is set to **"127.0.0.1"** and the Server Port is set to **"2400"**  
Click the **"Connect"** button

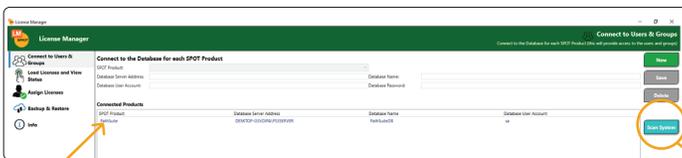


5 In the **"Connect to Users & Groups"** panel, press the **"Scan System"** button and confirm **"PathSuite"** is listed under **"Connected Products"**

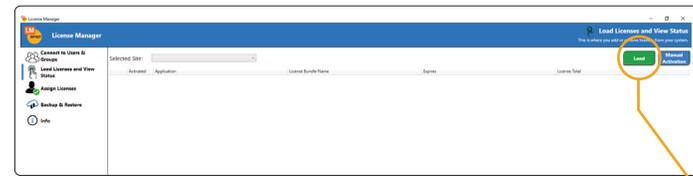

6 In the **"Load Licenses and View Status"** panel, click the **"Load"** button and select your **.lic** file



7 Wait for the license file to load and to automatically be activated (noted by a checked box in the **"Activated"** column)

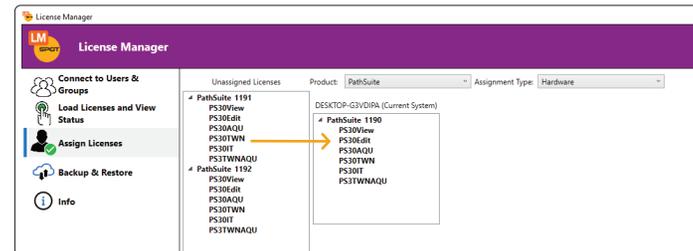

8 In the **"Assign Licenses"** panel, drag and drop a License from the **"Unassigned Licenses"** list to the **"Current System"**

9 Close License Manager